Operating Instructions

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Stirring Motors

Rührwerk

R 18, R 50, R 50D, R 80D-PC

Content

1	General Information	4
1.1	Scope of Delivery	4
2	User Instructions	5
2.1	Important Instructions for your safety	5
2.2	Danger symbols in this operating manual	5
2.2.1	Danger symbols:	5
2.2.2	Danger levels	5
3	General safety warnings and instructions	6
3.1	Warning signs on the device	7
4	Intended Use	7
4.1	General Description	7
5	Unpacking	8
6	Set-up	8
6.1	Mounting of Support Rod	8
6.2	Attach the Overhead Stirrer to a Stand	9
6.3	Attachment of Stirring Paddle	9
6.4	Assembling the Shaft Protector	10
7	Operation	10
7.1	Description of the front panels R 18 and R 50	11
7.2	Description of the front panels R 50D and R 80D-PC	11
7.2.1	Description of the rear R 80D-PC	12
7.3	Operation R 18, R 50 and R 50D	12
7.4	Operation R 80D-PC	12
7.5	Serial Port Operation (for R 80D-PC only)	13
7.5.1	Programming example:	14
8	Maintenance	15
9	Dismantling, Transport and Storage	16
9.1	Dismantling	16
9.2	Transport and Storage	16
10	Warranty and Liability	17
11	Technical Data	18
12	Disposal	19
13	Repairs	19
13.1	Repair Return Form	20
13	Konformitätserklärung / Declaration of conformity	38

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1 General Information

The overhead stirrers R 18, R 50, R 50D and R 80D-PC are designed in accordance with Safety Class 1, and built and tested in accordance with DIN EN 61010.

According to these regulations, the unit is designed to meet the requirements for safe and correct operations. To maintain the proper safety and operational functions of the instrument, the user should follow the instructions and safety guidelines in this manual.

1.1 Scope of Delivery

Inspect the content of the package for damage or missing parts:

	1 Overhead stirrer R 18	230 V (Part no.: 60214-0000)	with EU-plug
Or	1 Overhead stirrer R 18	230 V (Part no.: 60214-0003)	with UK-plug
Or	1 Overhead stirrer R 18	115 V (Part no.: 60214-0001)	with US-plug
Or	1 Overhead stirrer R 50	230 V (Part no.: 60215-0000)	with EU-plug
Or	1 Overhead stirrer R 50	230 V (Part no.: 60215-0003)	with UK-plug
Or	1 Overhead stirrer R 50	115 V (Part no.: 60215-0001)	with US-plug
Or	1 Overhead stirrer R 50D	230 V (Part no.: 60216-0000)	with EU-plug
Or	1 Overhead stirrer R 50D	230 V (Part no.: 60216-0003)	with UK-plug
Or	1 Overhead stirrer R 50D,	115 V (Part no.: 60216-0001)	with US-plug
Or	1 Overhead stirrer R 50D-SW,	230 V (Part no.: 60216-0013)	with UK-plug
Or	1 Overhead stirrer R 50D-SW,	115 V (Part no.: 60216-0011)	with US-plug
Or	1 Overhead stirrer R 80D-PC	230 V (Part no.: 60311-0000)	with EU-plug
Or	1 Overhead stirrer R 80D-PC	230 V (Part no.: 60311-0003)	with UK-plug
Or	1 Overhead stirrer R 80D-PC	115 V (Part no.: 60311-0001)	with US-plug
	1 Support Rod	(Part no.: 50215-0002)	
	2 Screws for Support Rod	(Part no.: 00215-0025)	
	1 Hexagon wrench key	(Part no.: 20285-0018)	

1 Instruction Manual

Accessories and Stirring Paddles (optional)

Guard for agitator shaft	(Part no.: 20618-0060)
Pedestal Stand incl. rod	(Part no.: 60491-0000)
Cross over clamp	(Part no.: 60492-0000)
Cross over clamp made of steel	(Part no.: 60492-0050)
Blade stirring paddle	(Part no.: 8B00940400)
Propeller stirring paddle	(Part no.: 8B00910500)

Optional Accessories for R 80D-PC

9 Pin SUB-D cable, 1.8 m long	(Part no.: 30275-0051)
USB-RS232 Adaptor	(Part no.: 30244-0001)

2 User Instructions

2.1 Important Instructions for your safety

- Every user must read and understand this manual completely before use. Only instructed users may operate the instrument. Failure to do so can result in serious injury or death.
- Follow general instructions for hazard prevention and general safety instructions, e.g. wear protection clothing, eye protection and gloves.
- This operating manual is part of the product. Thus, it must always be easily accessible.
- Enclose this operating manual when transferring the device to another place.
- If this manual is lost, please request another one. Please contact your dealer or

Ingenieurbüro CAT
M. Zipperer GmbH
Wettelbrunner Str. 6
D-79282 Ballrechten-Dottingen
Tel. ++49-7634-5056-800
info@cat-ing.de

2.2 Danger symbols in this operating manual

The safety instructions in this manual appear with the following danger symbols and danger levels:

2.2.1 Danger symbols:

	Hazard point		Electrical shock
	Risk of fire		Explosion
	Crushing		Bio hazard

2.2.2 Danger levels

 DANGER	Will lead to severe injuries or death
 WARNING	May lead to severe injuries or death
 CAUTION	May lead to light to moderate injuries
NOTICE	May lead to material damage

3 General safety warnings and instructions

	<p>⚠ DANGER Risk of explosion.</p> <ul style="list-style-type: none"> • Do not operate the device in the vicinity of highly flammable or explosive substances. The instrument is not explosion-proof. • Do not use this device for processing any substances which could generate an explosive atmosphere. • Do not use this device to process any explosive or highly reactive substances.
	<p>⚠ DANGER Electric shock as a result of penetration of liquid.</p> <ul style="list-style-type: none"> • Do not allow any liquids to penetrate the inside of the housing • Switch off the device and disconnect the power plug before starting cleaning or disinfection work. The On/Off Switch on the device does not disconnect the device from the power source. • Only plug the device back in if it is completely dry, both inside and outside.
	<p>⚠ WARNING Risk from incorrect supply voltage</p> <ul style="list-style-type: none"> • Only connect the device to an AC power source with a protective earth (PE). • Only connect the device to voltage sources which correspondent to the electrical requirements on the type label.
	<p>⚠ WARNING Electric shock due to damage to device or mains cable</p> <ul style="list-style-type: none"> • Only connect the device to the mains supply if the device and the mains cable are undamaged • Only use devices that have been properly installed or repaired. • In case of danger, disconnect the device from the mains supply by pulling the power plug from the mains socket or by using the isolating device intended for this purpose (e.g. emergency stop switch)
	<p>⚠ WARNING Lethal voltage inside the device</p> <ul style="list-style-type: none"> • Do not open the device. • Ensure that the housing is always closed and undamaged so that no parts inside the housing can be contacted by accident. • Do not allow any liquids to penetrate the inside of the housing. • Repairs are only to be carried out by trained service technicians.
	<p>⚠ WARNING Damages to health due to infectious liquids and pathogenic germs.</p> <ul style="list-style-type: none"> • When handling infectious liquids and pathogenic germs, observe the national regulations, the biological security level of your laboratory, the material safety data sheets and the manufacturer's application notes. • Wear personal protective equipment • For comprehensive regulations about handling germs or biological material of the risk group II or higher, please refer to the "Laboratory Biosafety Manual" in its respectively current valid version from the World Health Organisation
	<p>⚠ WARNING Risk of fire</p> <ul style="list-style-type: none"> • Do not use this device to process any highly flammable liquids
	<p>⚠ WARNING Damage to health due to contaminated device and accessories</p> <p>In the following cases, sample material can be released:</p> <ul style="list-style-type: none"> - improperly fixed agitator vessels - unstable vessels - damaged sealings - smashed glas vessels <ul style="list-style-type: none"> • Only mix in suitable vessels

	<ul style="list-style-type: none"> ● Observe the nationally prescribed safety environment when working with hazardous, toxic and pathogenic samples. Pay particular attention to personal protective equipment (gloves, clothing, goggles, etc.), extraction, and the safety class of the lab. ● Decontaminate the device and the accessories before storage and shipping.
	<p>CAUTION! Poor safety due to inadequate fixing of the unit</p> <ul style="list-style-type: none"> ● Ensure that the unit is firmly attached to a solid stand.
	<p>CAUTION! Poor safety due to incorrect accessories and spare parts.</p> <p>The use of accessories and spare parts other than recommended by Ingenieurbüro CAT, M. Zipperer GmbH may impair the safety, function and precision of the device. Ingenieurbüro CAT, M. Zipperer GmbH cannot be held liable or accept any liability for damage resulting from the use of incorrect or non-recommended accessories and spare parts, or from the improper use of such equipment.</p> <ul style="list-style-type: none"> ● Only use accessories and spare parts recommended by Ingenieurbüro CAT, M. Zipperer GmbH
	<p>CAUTION! Crush hazard due to moving parts</p> <ul style="list-style-type: none"> ● Always use a stirring shaft protector ● Do not replace any consumables as long as the device is running. ● Do not detach the stirring shaft protector as long as the device is running ● Only switch-on the device with properly mounted tools and completely closed stirring shaft protector.

3.1 Warning signs on the device

▲WARNING

This symbol indicates to read the instruction manual carefully prior to operation of the instrument. Please mark points which require special attention in your field of application so they are not overlooked. Disregarding of warnings may result in impairment of serviceability as well as impairment of the user.

4 Intended Use

The overhead stirrers R 18, R 50, R 50D and R 80D-PC are designed for use in chemical and biological laboratories of industrial enterprises, universities and pharmacies. They are suitable for stirring and mixing of liquids from aqueous to medium viscosity. For correct use the devices must be fixed to an appropriate stand.

To ensure maximum service life, observe the specified ambient conditions (temperature and humidity) and ensure that the instrument is not exposed to a corrosive atmosphere.

4.1 General Description

The laboratory overhead stirrers R 18, R 50 (without digital speed display), the R 50D (with digital speed display) and R 80D-PC are microprocessor controlled. They are particularly suitable for all applications where solutions of average viscosity have to be stirred. The use of the instrument is advantageous where the speed adjustment of stirring processes has to be reproducible - such as when test series have to be carried out one after the other under comparable conditions, and when parallel experiments with several stirrers must be made simultaneously.

The R 50, R 50D and R 80D-PC stirrers can also be used in the treatment of special solutions where higher viscosity values occur for short periods as the individual constituents are added.

The speed control is steplessly adjustable between 20 and 700 rpm for the R 80D-PC, between 50 and 1600 rpm for stirring motors R 50 and R 50D; for the R 18 the speed is adjustable between 110 and 2000 rpm. The electronic feedback control guarantees exact control and high precision so that fluctuations from the viscosity of the media being stirred will not occur.

5 Unpacking

Unpack the instrument carefully and check to see that it is not damaged. It is important that any damage incurred in transport be recognized at the time of unpacking. Notify your carrier or forwarding agent immediately in case of such damage.

▲WARNING

- Read this instruction manual carefully before operating the instrument. Should there be any additional questions, after reading these instructions, concerning the set-up, operation or warranty, please contact either your distributor, or the manufacturer.
- After reading and understanding the instruction manual you may now start operating the unit.
- Store the instruction manual in a place easily accessible to every user.

6 Set-up

▲DANGER ▲

- The units are not to be used in rooms with danger of explosion.
- Do not use this device to process any highly flammable liquids
- The unit is not to be used without supervision.


Please put the unit on to a fire-proof respectively non-combustible even surface.

6.1 Mounting of Support Rod

▲WARNING

The device operates with a high torque and vibrated when running. Insufficient mounted mechanical parts may detach themselves and endanger the health of the user.

- Ensure that all mechanical connections are firmly tightened.
- Ensure that the unit is firmly attached to a stand.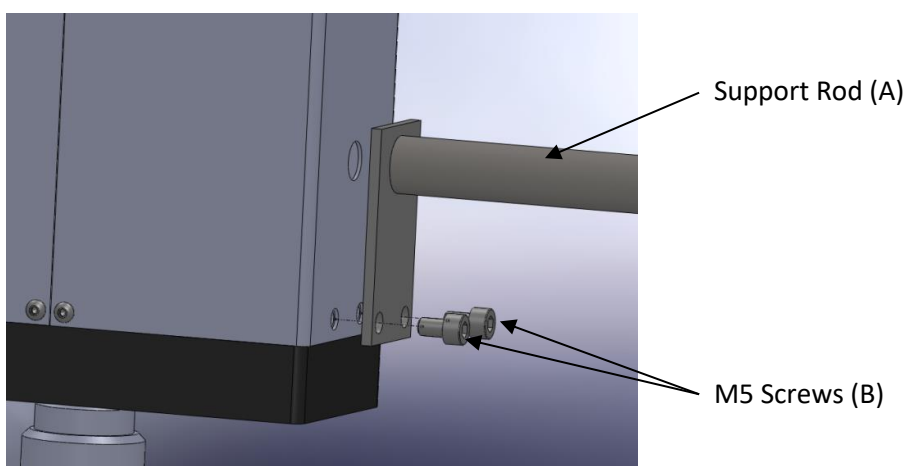

- Screw the support rod (A) to the rear panel of the overhead stirrer with the two M5 screws (B). Use a hexagon socket screw key with AF (across flats).
- Tighten the screws thoroughly.


6.2 Attach the Overhead Stirrer to a Stand

▲WARNING

Ensure that all mechanical connections are firmly tightened

Mount the cross over clamp (C) to the rod of the stand.

Insert the support rod (A) of the over head stirrer into the opening of the cross over clamp (C) and turn the knurled screw (D) until the drive unit is firmly attached to the stand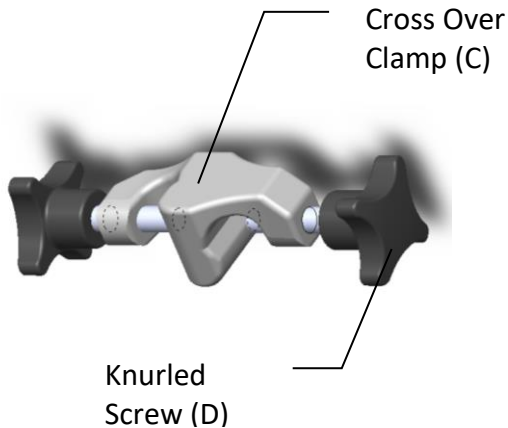


6.3 Attachment of Stirring Paddle

R 50, R 50D and R 80D-PC have a hollow shaft with a 10.5 mm bore. The chrome plated chuck avoids corrosion. The chuck takes any stirring rod up to 10 mm (stirring rod not included). The support rod at the back of the unit is used to attach the motor to a stand (stand not included).

R 18 does not have a hollow shaft but also takes any stirring rod up to 10 mm (stirring rod not included).

- Insert the stirring paddle into the chuck and firmly tighten it in the desired position by hand.
- The immersion depths of the stirring tool may be adjusted by loosening the chuck and axially moving the stirring tool.

6.4 Assembling the Shaft Protector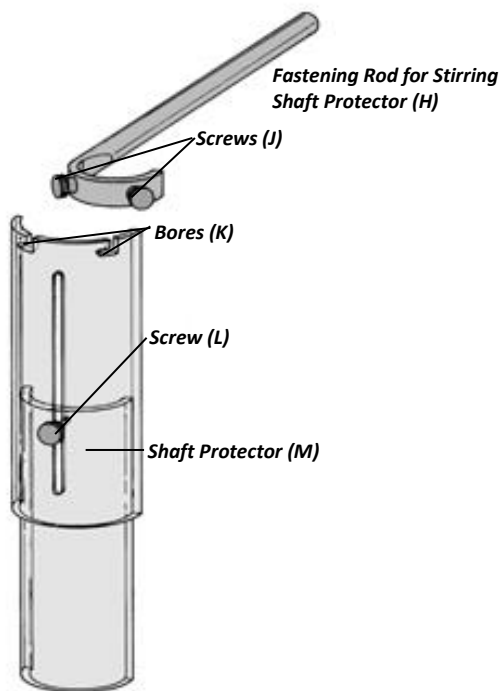

Use a stirring shaft protector to provide protection against injury when working with the unit.

Attach the fastening rod (H) of the shaft protector (M) in front of the chuck to the stand; e.g. use cross over clamp part no. 60492-0000) to fix the stirring shaft protector. Now bring the pockets (K) to the screws (J) and turn the stirring shaft protector to the left to the limit stop.

The screw (L) is used to adjust the length of the stirring shaft protector.

Check that the stirring shaft protector is held in position securely prior to each use and also at regular intervals. The position of the stirring shaft protector must only be adjusted when the unit is stationary and the power supply is disconnected.

7 Operation

⚠ DANGER

- The units are not to be used in rooms with danger of explosion.
- The unit is not to be used without supervision.

⚠ WARNING


- When connecting the instrument to an AC power outlet, ensure that your local supply voltage matches that indicated on the instrument's rating plate.
- Avoid splashing of liquids by using a suitable vessel, lower liquid level. Make sure that suitable protective clothing and eye-wear is used.
- This instruction sheet does not purport to address all of the safety problems which might result from the use of this instrument, chemicals, reagents, apparatus or equipment employed in any specific test or protocols. It is the responsibility of the user to consult their authorised safety advisors and establish appropriate health and safety practices and then determine the application of regulatory limitations prior to use.

Please put the unit on to a fire-proof respectively non-combustible even surface.


- Set the speed setting knob (7) to minimum.
- Press the power on/off button (9)
- Set the desired motor speed using the speed setting knob (7).

The digital value of the set motor speed can be read out in the top line of the display (1)

7.1 Description of the front panels R 18 and R 50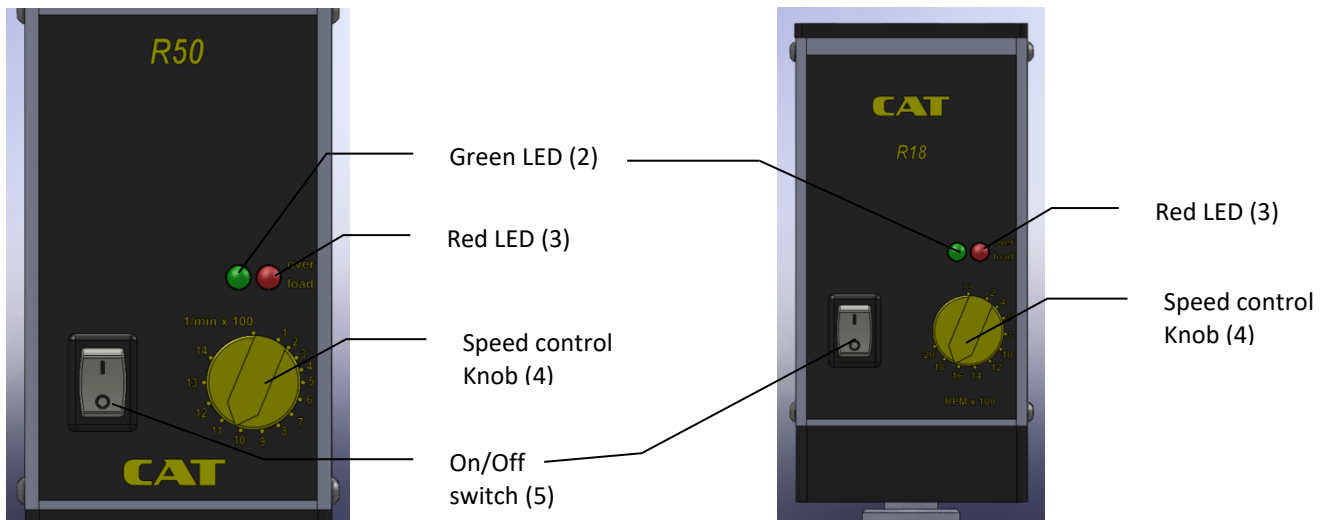


7.2 Description of the front panels R 50D and R 80D-PC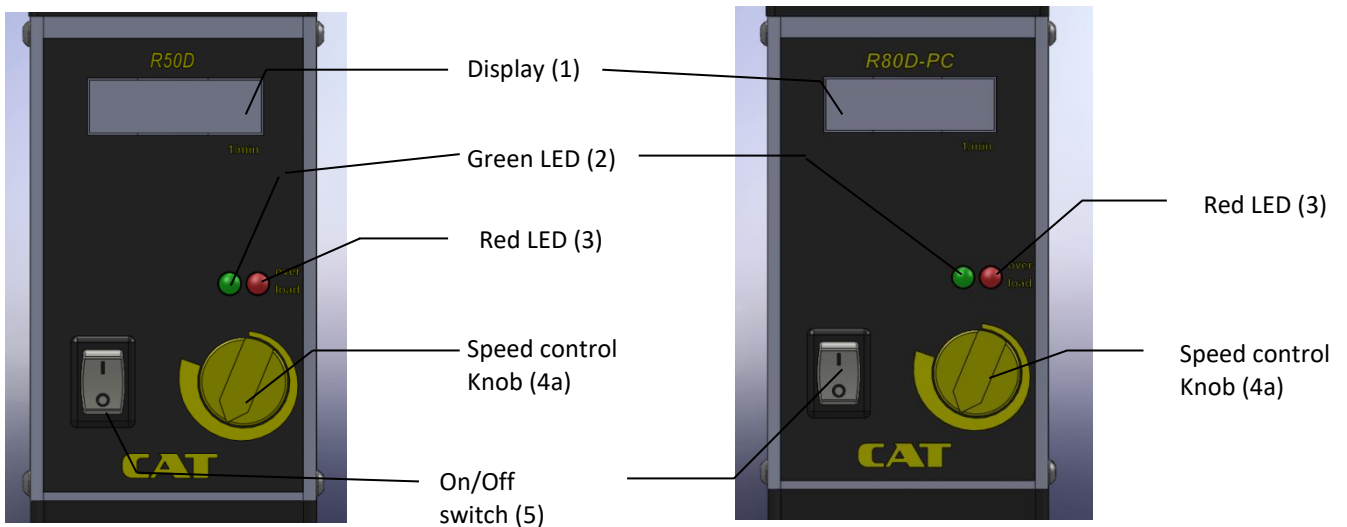

The R 50, R 50D and R 80D-PC stirrers have two independent safeguards against overload:

1. Should the allowable torque be exceeded, a red warning LED will show and the electronic system automatically reduces the power. As soon as the overload is either removed or overcome the red warning LED will go off. At continuous overload and the resulting reduced power to a speed $\text{rpm} < 20$ the motor will be switched off.
2. Should an overload occur from overheating during continual operations or too hot of an ambient, the red LED will show and the electronic will reduce the power intake. The motor will switch off if continuously overheated. After a cool down period and switching the unit off and on the motor will operate as normal.

7.2.1 Description of the rear R 80D-PC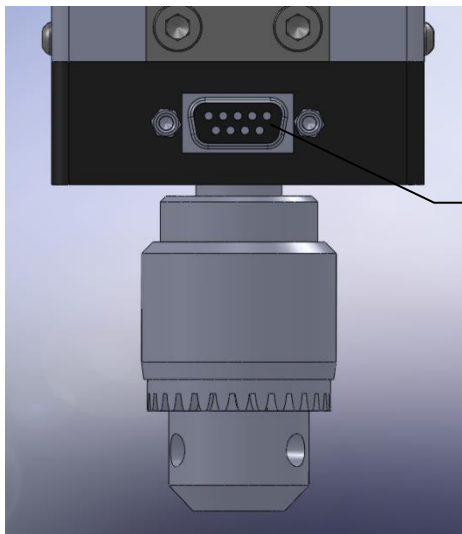

9-pin SUB- D
connection
RS232-interface (6)

7.3 Operation R 18, R 50 and R 50D

- Insert stirring rod into chuck and evenly tighten the chuck by hand.
- Insert plug into the mains outlet.
- Place stirring rod into the media to be stirred.
- Press switch to turn the device on.
- Adjust the speed by turning the speed regulator knob to the left. The speed for the R 18 and R 50 is reflected by the scale around the regulator knob. The speed for the R 50D is indicated on the digital display. Always start with the lowest speed.
- To turn off, reduce the speed to the lowest rate, turn the switch off, and then remove the stirring rod from the media.

7.4 Operation R 80D-PC


- Insert stirring rod into chuck and evenly tighten the chuck by hand.
- Insert plug into the mains outlet.
- Place stirring rod into the media to be stirred.
- Press switch to turn the device on.
- Adjust the speed by turning the speed regulator knob to the left. The speed for the R 80D-PC is indicated on the digital display. Always start with the lowest speed.
- Connect serial cable to the device. Set the regulator knob to the left end position (minimum RPM). Switch the device on to control the device over the serial port.
- To turn off, reduce the speed to the lowest rate, turn the switch off, and then remove the stirring rod from the media.

7.5 Serial Port Operation (for R 80D-PC only)

The following signals are led out on the 9-pin SUB-D connector on the rear side of the instrument:

- RS232 control signals

Pin Number	Description
1	n.c.
2	TxD (RS232 transmit data)
3	RxD (RS232 receive data)
4	n.c.
5	GND (signal ground)
6	n.c.
7	n.c.
8	n.c.
9	n.c.

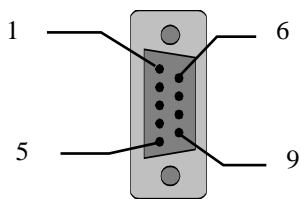*socket of the I/O connector (view on socket
– 9 pin Sub-D socket)*

Settings:

Baudrate: 1200
 Data bits: 8 bit
 Parity: none
 Stop bits: 1

Commands:

Query of version 0,RTY,1
 Change operation mode 0,WSM,1 (1 – PC, 0 – manual)
 Set RPM: 0,WSE,xxx (“xxx” = 20 to 700)
 Read RPM: 0,RSE,1
 Write device address: 0,WSA,x (“x” = 1 to 8)

7.5.1 Programming example:

Example 1: set RPM

1,WSM,1 ;PC mode
1,WSE,100 ;programs speed, (device address "1")

The device replies with the following code string:

1,HS,OK ;command has been accepted and executed

Example 2: Read RPM

1,WSM,1 ;PC mode
1,RSE,1 ;Read RPM of device 1

The device replies with the following code string:

1,HS,OK,100 ;command has been accepted and executed; current RPM is 100

8 Maintenance

WARNING

Do not open the instrument. Repairs are only to be carried out by trained service technicians.

Apart from the motor carbon brushes the instruments are maintenance-free, except from normal wear and the statistical failure rate of the components. Should the instrument require repairs, it is recommended that they be done at the factory. Please send any defective instrument either to the Ingenieurbüro CAT M. Zipperer GmbH factory or your agent.

WARNING

Unplug the AC power cord before opening the housing.

The carbon brushes in the motor are subject to natural wear. To change the carbon brushes please carry out the following steps.

1. Remove the black cover at the side of the motor.
2. Bent up the metal clip and take out the spring.
3. Use a small screwdriver to remove the clip from the carbon and then take out the carbon.
4. Replace the carbon. Attach the spring. Bent aback the clip. Attach the black cover.

Should the tooth belt have to be changed, then:

NOTICE

The following information does not apply for overhead stirring motor R 80D-PC!!!

1. Remove the hard plastic undercover.
2. Pull the tooth belt from the large cogwheel.
3. Put the new tooth belt on the small cogwheel and then onto the large cogwheel.
4. Now turn lightly the large cogwheel until the belt is in the proper position.

All bearings of the stirring motor are sufficiently lubricated so that with careful handling no maintenance is necessary. Should problems occur, please send the instrument to Ingenieurbüro CAT M. Zipperer GmbH or your agent.

9 Dismantling, Transport and Storage

9.1 Dismantling

1. Switch the unit off.
 2. Disconnect the unit from the mains.
-
- Remove any glass beakers and any other equipment around the overhead stirrer.
 - Remove the stirring tool by loosening the chuck
 - Loosen the cross over clamp and remove the clamp from the support rod
 - Unscrew the support rod.
 - Now you may remove the instrument from the working area.

9.2 Transport and Storage

Prior to transport:

Switch the instrument off and unplug from power supply.

Place the instrument and its parts in its original packaging or another suitable container to protect it during transport. Close the packaging with adhesive tape.

Store the instrument in a dry environment. Please observe the specified conditions of the ambient (temperature and humidity).

Do not subject the instrument to mechanical shocks or vibration during transporting it.

In case you do not use the original packaging please mark the box with the following notes:

- Glass symbol (handle with care, fragile)
- Umbrella (keep dry)
- Content (list of content)

10 Warranty and Liability

The manufacturer agrees to correct for the original user of this product, either by repair, or at the manufacturer's election, by replacement, any defects in material or workmanship which develop within 24 months after delivery of this product to the original user. In the event of replacement, the replacement unit will be warranted for the remainder of the original twenty-four (24) months period of ninety (90) days, whichever is longer.

If this product should require service, contact your distributor or manufacturer for necessary instructions.

This warranty shall not apply if the defect or malfunction was caused by accident, neglect, unreasonable use, improper service, or other causes not arising out of defects in material or workmanship.

There are no warranties, expressed or implied, including, but not limited to, those of merchantability or fitness for a particular purpose, which extended beyond the description and period set forth herein.

The manufacturer's sole obligation under this warranty is limited to the repair or replacement of a defective product and the manufacturer shall not, in any event, be liable for any incidental or consequential damages of any kind resulting from use or possession of the product.

11 Technical Data

R 80D-PC	
Motor	Permanently generated direct –current motor
Speed Range	20 – 700 1/min
Nominal Rating	120 Watt input Short Term Overload 30%
Torque over entire speed range	80 Ncm
Capacity of Chuck	0.5 – 10.5 mm
Viscosity of Medium	Average values (water to wallpaper paste)
Mains Voltage	200 -250 V, 50/60 Hz or 115 V +/- 10%, 50/60 Hz
Permissible period of operation	Continuous operation may require service intervals
Permissible ambient temperature	+ 5°C to 40°C
Diameter of support rod	16 mm
Length of support rod	196 mm
Weight	2.5 kg
Dimensions	65 x 200 (incl. 55 mm chuck) x 120 mm (w x h x d)

R 50 and R 50D	
Motor	Permanently generated direct –current motor
Speed Range	50 – 1600 1/min
Nominal Rating	120 Watt input Short Term Overload 30%
Torque over entire speed range	55 Ncm
Capacity of Chuck	0.5 – 10.5 mm
Viscosity of Medium	Average values (water to wallpaper paste)
Mains Voltage	200 -250 V, 50/60 Hz or 115 V +/- 10%, 50/60 Hz
Permissible period of operation	Continuous operation may require service intervals
Permissible ambient temperature	+ 5°C to 40°C
Diameter of support rod	16 mm
Length of support rod	196 mm
Weight	2.3 kg
Dimensions	65 x 200 (incl. 55 mm chuck) x 120 mm (w x h x d)

R 18	
Motor	Permanently generated direct –current motor
Speed Range	100 – 2000 1/min
Nominal Rating	120 Watt input Short Term Overload 30%
Torque over entire speed range	18 Ncm
Capacity of Chuck	0.5 – 10.5 mm
Viscosity of Medium	Aqueous fluids
Mains Voltage	200 -250 V, 50/60 Hz or 115 V +/- 10%, 50/60 Hz
Permissible period of operation	Continuous operation may require service intervals
Permissible ambient temperature	+ 5°C to 40°C
Diameter of support rod	16 mm
Length of support rod	196 mm
Weight	2.2 kg
Dimensions	65 x 200 (incl. 55 mm chuck) x 120 mm (w x h x d)

12 Disposal

Please dispose of used instruments and defective components at your local recycling collection point. Prior to disposal, sort according to materials: metal, glass, plastic, etc. Also be sure to dispose of the packing material in an environmental-friendly manner.

13 Repairs

When returning instruments for repair that have come into contact with hazardous substances, please:

Fill in attached "Repair Return Form"

Provide precise information on the relevant medium

Take protective measures to ensure the safety of our receiving and maintenance personnel

Mark the package as appropriate for hazardous materials.

13.1 Repair Return Form

(Please copy and enclose the return)

Important:

1. Calibration of the device is done automatically.
2. You will receive an estimate for the expected repair costs.
3. Repairs will be done after release.

CONTACT/USER INFORMATION

Contact:	Phone No.
Fax No.	Email:
Billing:	Shipping:
Company	Company
Address	Address

DEVICE INFORMATION

Model	Serial No.
Please describe all problems/malfunctions	

Operating Conditions (please fill in if applicable)		
Ambient Temperature	Humidity	Speed
Load	Volume	Viscosity
Temperature in °C	Sample Temperature	Operating Time
Sample Description*		

*NOTE: If the device was exposed to hazardous material, it must be decontaminated BEFORE returning it to Ingenieurbüro CAT M. Zipperer GmbH and an MSDS for hazardous material must be included with the device.

RETURN SHIPPING

UPS <input type="checkbox"/>	Air Parcel Post <input type="checkbox"/>	Collect** <input type="checkbox"/>	Other** <input type="checkbox"/>
------------------------------	--	------------------------------------	----------------------------------

**Your account number is required for UPS collect respectively the address and contact of your preferred forwarder if you choose any other transport means

PACKAGING INSTRUCTIONS TO RETURN A DEVICE FOR REPAIR

- ✓ Remove all accessories (e.g. homogenizer tools, stirring paddles) from the device
- ✓ Clean excess testing material off the device/accessory
- ✓ Include MSDS for all hazardous materials used with this device.
- ✓ Pack the device in its original box. If the box is not available, take care to wrap the device and accessories with enough material to support them.
- ✓ Do NOT send pedestal stand unless there is a problem with the upright rod, clamp or base. If there is a problem with the stand remove the upright rod from the base and individually wrap each item to avoid contact with the device. (Applicable for overhead stirrers and homogenisers)
- ✓ Pack the device and related items in a strong box for shipping. Mark the outside of the box with handling instructions

Example: "Handle with care" or "Fragile-Delicate Instrument" and send to:

Ingenieurbüro CAT, M. Zipperer GmbH, Service Department
Wettelbrunner Str. 6
D-79282 Ballrechten-Dottingen

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We declare under our sole responsibility that this product (see table) corresponds to the regulations and conforms with the standards or standardized documents:

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The agreement with further valid guidelines/regulations following for the product is explained:

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Applied one harmonized standards in particular:

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In the case of a modification of the unit which has not been agreed on with us, this declaration becomes null and void and the warranty expires.

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