CAT

**Ingenieurbüro CAT
M. Zipperer GmbH**

Operation Instructions

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Stirring Motor

Rührmotor

R 16

Content

1	User Instructions	4
1.1	Important Instructions for your safety	4
1.2	Danger symbols in this operating manual.....	4
1.2.1	Danger symbols:.....	4
1.2.2	Danger levels.....	5
1.3	Warning signs on the device	5
2	General safety warnings and instructions	5
3	General Information	7
4	Unpacking.....	7
4.1	Scope of delivery	7
5	Intended Use.....	8
6	Operation	8
6.1	Description of front panel	8
6.1.1	Operating control.....	8
6.2	Assembling of support rod on the device	8
6.3	Mounting the stirring tool.....	9
6.4	Electrical Connection.....	9
7	Maintenance	9
8	Dismantling, Transport and Storage	10
8.1	Dismantling.....	10
8.2	Transport and Storage.....	10
9	Warranty	10
10	Technical Data.....	11
11	Disposal	11
11.1	Repair Return Form	12
22	Declaration of conformity.....	22

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1 User Instructions

1.1 Important Instructions for your safety

- Every user must read and understand this manual completely before use. Failure to do so can result in serious injury or death.
- Comply with all safety and accident-prevention regulations applicable to laboratory work.
- Follow general instructions for hazard prevention and general safety instructions, e.g. wear protection clothing, eye protection and gloves.
- This operating manual is part of the product. Thus, it must always be easily accessible.
- Enclose this operating manual when transferring the device to another place.
- If this manual is lost, please request another one. Please contact your dealer or

Ingenieurbüro CAT
M. Zipperer GmbH
Wettelbrunner Str. 6
D-79282 Ballrechten-Dottingen
Tel.: ++49-(0)7634-5056-800
Fax: ++49-(0)7634-5056-801
www.cat-ing.de
info@cat-ing.de

1.2 Danger symbols in this operating manual

The safety instructions in this manual appear with the following danger symbols and danger levels:

1.2.1 Danger symbols:

 Hazard point	 Explosion
 Risk of fire	 Chemical hazard
 Crushing	 Bio hazard

1.2.2 Danger levels

	Will lead to severe injuries or death
	May lead to severe injuries or death
	May lead to light to moderate injuries
	May lead to material damage

1.3 Warning signs on the device

WARNING

This symbol indicates that it is imperative to read and understand the instruction manual prior to operating the instrument. Please highlight points which require special attention in your field of application so they are not overlooked. Disregard of warnings may result in impairment of serviceability as well as in physical harm to the user.

2 General safety warnings and instructions

	<p> DANGER Risk of explosion.</p> <ul style="list-style-type: none"> • Do not operate the device in the vicinity of highly flammable or explosive substances. The instrument is not explosion-proof. • Do not use this device for processing any substances which could generate an explosive atmosphere. • Do not use this device to process any explosive or highly reactive substances. • Do not use flammable or explosive substances near the instrument.
	<p> WARNING Damage to health due to corrosive or aggressive chemicals</p> <ul style="list-style-type: none"> • Observe all markings on the reagent bottles. • Always check the instrument for leaks and air bubbles. Special attention should be directed to determine that all push-ons, threaded connections and suction tubes are firmly in place before beginning operation. Leaking solutions may endanger persons and materials • Use proper connecting vessels, protective clothing and gloves. • Avoid splashes • When dispensing, maintain a physical distance between the instrument and the body. • Dangerous and fuming chemicals must be dispensed in a fume hood. • Only employ the instrument for the purpose intended by the manufacturer, and particularly within the resistance limits of the instrument. If in doubt, contact your supplier, or the manufacturer's factory representative at the phone number shown at the front page of this operating instruction. • Always use the instrument in such a manner that neither the operator, nor any other person is endangered.
	<p>Damages to health due to infectious liquids and pathogenic germs.</p>

	<p>▲WARNING ● When handling infectious liquids and pathogenic germs, observe the national regulations, the biological security level of your laboratory, the material safety data sheets and the manufacturer’s application notes.</p> <ul style="list-style-type: none"> ● Wear personal protective equipment ● For comprehensive regulations about handling germs or biological material of the risk group II or higher, please refer to the “Laboratory Biosafety Manual” in its respectively current valid version from the World Health Organisation
	<p>▲WARNING Damage to health due to contaminated device and accessories</p> <p>In the following cases, sample material can be released:</p> <ul style="list-style-type: none"> - improperly sealed tubes - unstable tubes - high vapour pressure of the content so that the seal of the tubes can spring open - damaged sealing - smashed glass tubes <ul style="list-style-type: none"> ● Observe the nationally prescribed safety environment when working with hazardous, toxic and pathogenic samples. Pay particular attention to personal protective equipment (gloves, clothing, goggles, etc.), extraction, and the safety class of the lab. ● Decontaminate the device and the accessories before storage and shipping.
	<p>▲WARNING Risk of fire</p> <ul style="list-style-type: none"> ● Do not use this device to process any highly flammable liquids
	<p>▲CAUTION Poor safety due to incorrect mounted stirring paddle.</p> <p>For safe operation, a minimum clamping length of 80 mm is required.</p>
	<p>▲CAUTION Poor safety due to inadequate fixing of the unit</p> <ul style="list-style-type: none"> ● Ensure that the unit is firmly attached to a solid stand.
	<p>▲CAUTION Poor safety due to incorrect accessories and spare parts.</p> <p>The use of accessories and spare parts other than recommended by Ingenieurbüro CAT, M. Zipperer GmbH may impair the safety, function and precision of the device. Ingenieurbüro CAT, M. Zipperer GmbH cannot be held liable or accept any liability for damage resulting from the use of incorrect or non-recommended accessories and spare parts, or from the improper use of such equipment.</p> <ul style="list-style-type: none"> ● Only use accessories and spare parts recommended by Ingenieurbüro CAT, M. Zipperer GmbH
	<p>▲CAUTION Crush hazard due to moving parts</p> <ul style="list-style-type: none"> ● Do not replace any consumables as long as the device is running. ● Do not open the coverage as long as the device is running

3 General Information

The overhead stirrer R 16 is designed in accordance with Safety Class 1, and built and tested in accordance with DIN EN 61010.

According to these regulations, the unit is designed to meet the requirements for safe and correct operations. To maintain the proper safety and operational functions of the instrument, the user should follow the instructions and safety guidelines in this manual.

	<p>▲WARNING <i>Read this instruction manual carefully before operating the instrument. Should there be any additional questions, after reading these instructions, concerning the set-up, operation or warranty, please contact either your distributor, or the manufacturer at the following address:</i></p>
---	---

Ingenieurbüro CAT
M. Zipperer GmbH
Wettelbrunner Str. 6
D-79282 Ballrechten-Dottingen
Tel.: ++49-(0)7634-5056-800
Fax: ++49-(0)7634-5056-801
www.cat-ing.de
info@cat-ing.de

4 Unpacking

Unpack the instrument carefully and check to see that it is not damaged. It is important that any damage incurred in transport be recognised at the time of unpacking. Notify your carrier or forwarding agent immediately in case of such damage.

4.1 Scope of delivery

1 Stirring Motor R 16	part no.: 60213-0050
1 Support Rod	part no.: 10402-0064
1 Operating manual	
1 Power Supply	
consisting of:	
Switching Power Supply 24V DC	part no.: 30213-0054
Set of input plugs	part no.: 30213-0057

	<p>▲WARNING For your safety: Follow the instructions supplied with the power supply.</p>
---	--

Optional Accessories:

Pedestal stand,	part no.: 60491-0000
Cross over clamp,	part no.: 60492-0000 (aluminium), 60492-0050 (CRES)
Propeller Stirrer Ø6x300mm 2 blades 50 mm	part no.: 8B00970000
Propeller Stirrer Ø6x300mm 3 blades 50 mm	part no.: 8B00910000
Propeller Stirrer Ø6x300mm 4 blades 50 mm	part no.: 8B00900000
Stirring shaft Ø6x300mm, shaft end with M6, 10 mm long, without impeller	part no.: 10618-0001

5 Intended Use

The overhead stirrer R 16 is designed for use in chemical and biological laboratories of industrial enterprises, universities and pharmacies. They are suitable for stirring and mixing of liquids from aqueous to medium viscosity. For correct use, the devices must be fixed to an appropriate stand.

To ensure maximum service life, observe the specified ambient conditions (temperature and humidity) and ensure that the instrument is not exposed to a corrosive atmosphere.

6 Operation

	⚠ DANGER <i>The unit may not be operated in explosion-prone areas.</i>
	⚠ CAUTION <i>Pull main plug as a protection against unauthorised or accidental operation whenever the work is interrupted for extended periods.</i>

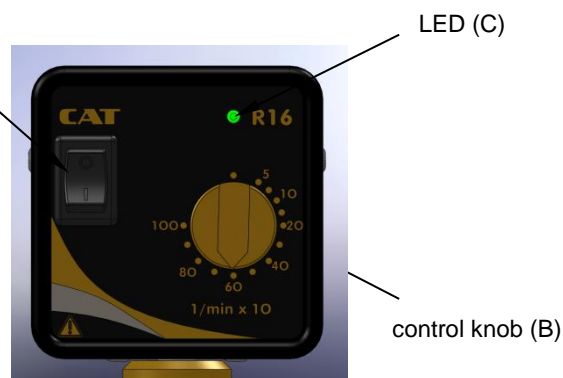
6.1 Description of front panel

6.1.1 Operating control

The following operation controls are available

- rocker switch (A) – switch on/off
- the control knob (B) is also for speed control
- LED (C) as status index:
 - green: set speed is reached
 - orange: set speed is not reached, load is high
 - orange flashing: overload – motor is switched off – remove load – restart via switching off/on

Power switch (A)

LED (C)

control knob (B)

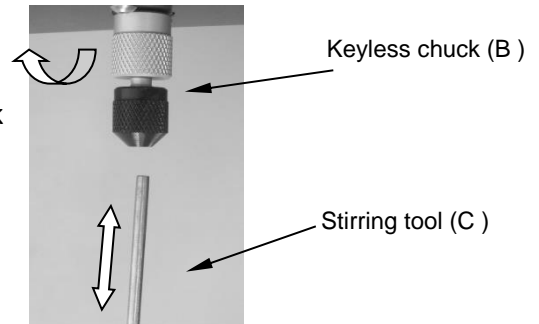
6.2 Assembling of support rod on the device

- Screw the support rod to the rear panel of the housing (clockwise).
- Tighten the support rod strongly by using a suitable 6 mm tool.
Max. thread length 7mm!

6.3 Mounting the stirring tool

The stirrer R 16 is a convenient, robust overhead stirrer for use with $\varnothing 6$ mm stirring rods.

1. Unfasten the keyless chuck (B)
2. Move the piston of the stirring tool (C) into the chuck
3. Torque down the chuck (B) when the piston is in the intended position

	CAUTION Poor safety due to incorrect mounted stirring paddle. For safe operation, a minimum clamping length of 80 mm is required.
--	--

6.4 Electrical Connection

	WARNING <i>The unit may not be operated in explosion-prone areas.</i>
	WARNING <i>Always start with the lowest speed. Increase the speed until the speed slowly until the desired speed setting has been reached.</i>

1. Insert plug into the mains outlet.
2. Place stirring rod into the media to be stirred.
3. Turn the device on via rocker switch (A)

The drive motor is designed for continuous operation.
It has an electronically controlled speed from 50 - 1000 rpm.

7 Maintenance

NOTICE *Do not open the instrument. Repairs are only to be carried out by trained service technicians.*

This instrument is guaranteed for a period of 24 months in accordance with our terms of warranty. Please see warranty for details. In case of malfunction, do not attempt to repair the unit. Any attempt by the user to repair the unit will cancel the warranty. Contact the distributor.

Due to safety and the warranty, the Stirrer R 16 should only be opened and repaired by authorised service personnel. Any work on the electronics of the unit should only be done by knowledgeable and trained personnel. Any attempt by the user to repair the unit will cancel the warranty.

All bearings of the stirring motor are sufficiently lubricated so that with careful handling no maintenance is necessary. Should problems occur, please send the instrument to Ingenieurbüro CAT, M. Zipperer GmbH or your agent.

	WARNING <i>Before opening the device unplug the mains plug.</i>
--	--

8 Dismantling, Transport and Storage

8.1 Dismantling

1. Switch the unit off.
2. Disconnect the unit from the mains.
3. Remove any glass beakers and any other equipment around the overhead stirrer.
4. Remove the stirring tool by loosening the chuck
5. Loosen the cross over clamp and remove the clamp from the support rod
6. Unscrew the support rod.
7. Now you may remove the instrument from the working area.

8.2 Transport and Storage

Prior to transport:

Switch the instrument off and unplug the power supply.

- Place the instrument and its parts in its original packaging or another suitable container to protect it during transport. Close the packaging with adhesive tape.
- Store the instrument in a dry environment. Please observe the specified conditions of the ambient (temperature and humidity).
- Do not subject the instrument to mechanical shocks or vibration during transporting it.
- In case you do not use the original packaging please mark the box with the following notes:
 - Glass symbol (handle with care, fragile)
 - Umbrella (keep dry)
 - Content (list of content)

9 Warranty

The manufacturer agrees to correct for the original user of this product, either by repair, or at the manufacturer's election, by replacement, any defects in material or workmanship which develop within 24 months after delivery of this product to the original user. In the event of replacement, the replacement unit will be warranted for the remainder of the original 24 months period of ninety (90) days, whichever is longer.

If this product should require service, contact your distributor or manufacturer for necessary instructions.

This warranty shall not apply if the defect or malfunction was caused by accident, neglect, unreasonable use, improper service, or other causes not arising out of defects in material or workmanship.

There are no warranties, expressed or implied, including, but not limited to, those of merchantability or fitness for a particular purpose, which extended beyond the description and period set forth herein.

The manufacturer's sole obligation under this warranty is limited to the repair or replacement of a defective product and the manufacturer shall not, in any event, be liable for any incidental or consequential damages of any kind resulting from use or possession of the product.

10 Technical Data

Part No.	60213-50
Power supply	Switching Power Supply, Input 100-240V, Output 24V , 25W
Stirring Volume (H₂O)	10 l
Motor	brushless DC Motor, electronically controlled speed
Speed Range	50 - 900 1/min
Torque	4 Ncm
Weight incl. Power supply	approx. 0.7 kg.
Mounting support	188 x Ø12 mm
Chuck	Ø 6 mm
Dimensions	67 x 67 x 121 mm plus chuck 38mm x Ø18mm

Conditions of surroundings:

Admissible temp. of surrounding	+ 5°C up to + 40°C max. 80 % relative humidity
Admissible operation	continuous operation

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▲WARNING *When returning instruments for repair that have come into contact with hazardous substances, please:
Fill in attached "Repair Return Form"
Provide precise information on the relevant medium
Take protective measures to ensure the safety of our receiving and maintenance personnel
Mark the package as appropriate for hazardous materials.*

11.1 Repair Return Form

Important: <ol style="list-style-type: none"> 1. Calibration of the device is done automatically. 2. You will receive an estimate for the expected repair costs. 3. Repairs will be done after release.

CONTACT/USER INFORMATION	
Contact:	Phone No.
Fax No.	Email:
Billing:	Shipping:
Company	Company
Address	Address

DEVICE INFORMATION	
Model	Serial No.
Please describe all problems/malfunctions	

Operating Conditions (please fill in if applicable)		
Ambient Temp.	Humidity	Speed
Load	Volume	Viscosity
Temperature in °C	Sample Temperature	Operating Time
Sample Description*		

*NOTE: If the device was exposed to hazardous material, it must be decontaminated BEFORE returning it to Ingenieurbüro CAT, M. Zipperer GmbH and an MSDS for hazardous material must be included with the device.

RETURN SHIPPING			
UPS <input type="checkbox"/>	Air Parcel Post <input type="checkbox"/>	Collect** <input type="checkbox"/>	Other** <input type="checkbox"/>

**Your account number is required for UPS collect respectively the address and contact of your preferred forwarder if you choose any other transport means

PACKAGING INSTRUCTIONS TO RETURN A DEVICE FOR REPAIR

- ✓ Remove all accessories (e.g. homogenizer tools, stirring paddles) from the device
- ✓ Clean excess testing material off the device/accessory
- ✓ Include MSDS for all hazardous materials used with this device.
- ✓ Pack the device in its original box. If the box is not available, take care to wrap the device and accessories with enough material to support them.
- ✓ Do NOT send pedestal stand unless there is a problem with the upright rod, clamp or base. If there is a problem with the stand remove the upright rod from the base and individually wrap each item to avoid contact with the device. (Applicable for overhead stirrers and homogenisers)
- ✓ Pack the device and related items in a strong box for shipping. Mark the outside of the box with handling instructions
Example: "Handle with care" or "Fragile-Delicate Instrument" and send to:

Ingenieurbüro CAT, M. Zipperer GmbH, Wettelbrunner Str. 6, D-79282 Ballrechten-Dottingen

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We declare under our sole responsibility that this product (see table) corresponds to the regulations and conforms with the standards or standardized documents:

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The agreement with further valid guidelines/regulations following for the product is explained:

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Applied one harmonized standards in particular:

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In the case of a modification of the unit which has not been agreed on with us, this declaration becomes null and void and the warranty expires.

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User Instructions Power supply GS-Series

Foreword

This AC/DC adapter is designed as external power supply for electrical equipment. Please read and follow the instructions before first use.

Installation

Before first use please check the mains voltage to match the input voltage specified on the power supply. Then check the output voltage, its polarity, the output current and power to match the requirements of the equipment to be connected with.

The connection of the plug must be stable; twisting or bending of the cable should be avoided.

Provide adequate ventilation during operation. The adapter should not be covered or used near the heating sources and may not be exposed to direct sunlight.

If your device is not used for a long time we recommend to disconnect the adaptor from main voltage in order to avoid any damage due to overvoltage for example (spikes, lightning).

Safety

Do not use the adapter if it is open or its housing is damaged; furthermore is prohibited to continue using the adapter in case of main connection damage - life danger due to potentially electric shock.

Repairs may be performed only by an authorized technician; any unauthorized modifications or repairs to the unit will void the warranty.

Make sure the DC-connector on secondary side of the adapter is suitable for your device - any forced connection of a non-matching DC connector may damage the device.

Protect the adapter from water and moisture.

Maintenance

For cleaning you may not use cleaning agents. Wipe with a dry cloth. Disconnect the adapter from the mains before cleaning.

Disposal

Electronic devices are hazardous waste and do not belong to the normal house waste. If the device reached end of its life dispose it according to the applicable legal provisions of the municipal collection points. Disposing in the domestic waste is not permitted.

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MEAN WELL ENTERPRISE Co., Ltd.
Schukat electronic Vertriebs GmbH